



# Quickstart Installer Guide

## Complete Stand-alone Network Connected System Kit


Remove the touch screen computer, keyboard, mouse, and power supply from the box. The optional Barcode Scanner is in a separate box.

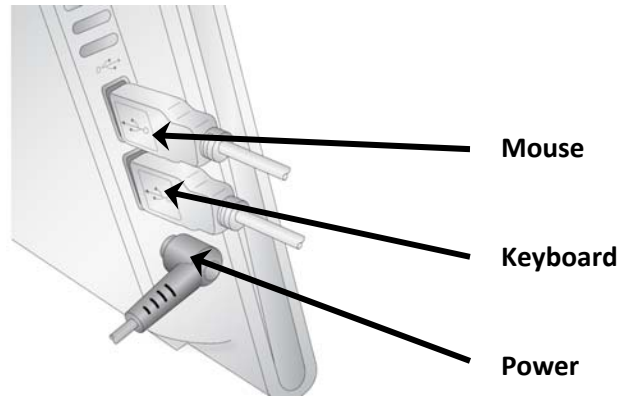

1. Connect the keyboard, mouse and power connector to the left side of the touch screen computer.

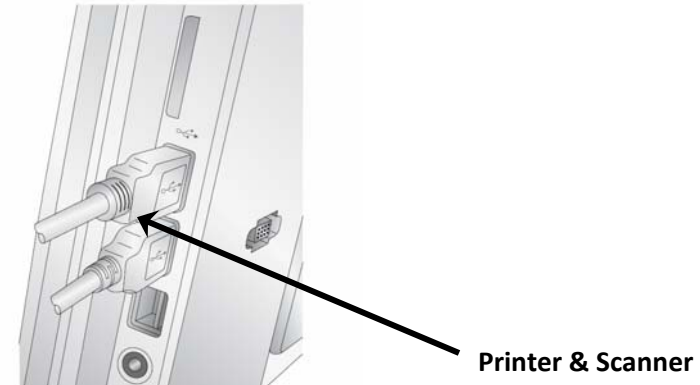

2. Connect the USB cable for the Slide Printer to the right side of the touch screen computer and the back of the printer. Connect the Scanner (optional) to the right side of the computer.

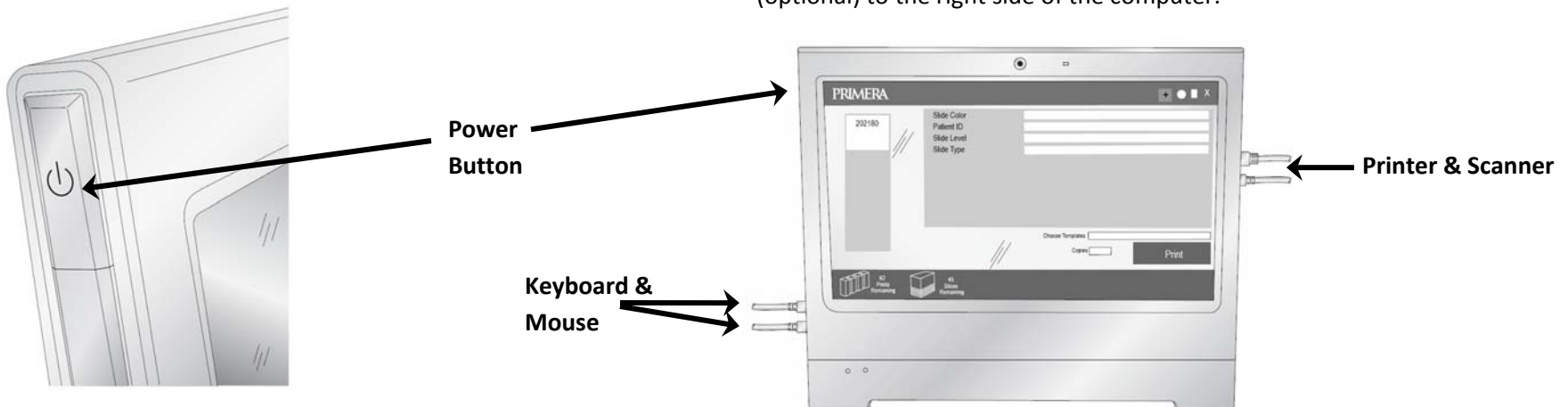

3. Press the power button on the top left side of the touch screen computer.

4. PTLab™ will automatically start. (If it does not start or if it is accidentally closed you can double click the PTLab Icon on the desktop to restart it.)